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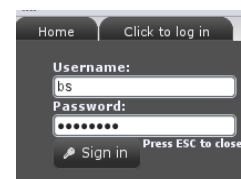
14. Klik på "Send".

The screenshot shows the first step of a three-step process. The top navigation bar has three tabs: '1 Login' (selected), '2 Attachments', and '3 Finalize'. The main content area is titled 'Login & select slide box...' and contains two checked status items: 'You are logged in as bs' and 'Your current slidebox is auhisto'. At the bottom, there are three buttons: 'Previous', 'Next' (highlighted), and 'Close'.

The screenshot shows the second step of the process. The top navigation bar has three tabs: '1 Login', '2 Attachments' (selected), and '3 Finalize'. The main content area is titled 'Step 2 Content' and includes a 'Click to refresh...' link. Below this, there are two columns of text. The left column lists '1 Lymph node', 'Retikulære fibre', 'fedtvæv', 'Mitosis', and 'Reticulum celle'. The right column lists '10 Bone marrow' and 'mitosis'. A vertical scrollbar is visible on the right side of the content area. At the bottom, there are three buttons: 'Previous', 'Next' (highlighted), and 'Close'.

The screenshot shows the third step of the process. The top navigation bar has three tabs: '1 Login', '2 Attachments', and '3 Finalize' (selected). The main content area is titled 'Step 3 Content' and contains a 'Recipient's email:' label. Below this are two input fields: the first contains 'bsander@ki.au.dk;navn@eks.dk' and the second contains 'HistoViewer Test'. A large text area for the message body is located below the input fields. At the bottom right of the text area is a 'Send' button with an envelope icon. At the bottom of the entire window, there are three buttons: 'Previous', 'Next', and 'Close' (highlighted).